Easy UV

Application Software

For UV/VIS Spectrophotometers

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Tel: +39 059 653 274

Please contact your local distributor

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ENGLISH

1. Feature overview

Easy UV is the application software compatibles with Microsoft Windows® operating systems for the remote control of the UV/VIS ONDA Spectrophotometers TOUCH Series. The main operating functions of the Easy UV software are summarized below.

Quantitative Analysis

Measuring the concentration of adequately prepared sample through a calibration curve. There are two ways to construct a calibration curve:

- 1) measurements of Standard solutions with a known concentration (up to 20 including repeated tests)
- 2) direct insertion of the coefficients (factors)

Both modes above include three types of regression:

- 1) Linear
- 2) Linear through origin
- 3) Quadratic

Kinetics Analysis

Measuring Absorbance or Transmittance of a sample over time at regular intervals of time (selectable between 0.5 / 1/2/5/10/30/60 seconds).

The Absorbance/time graph and the data table are displayed on the screen.

Spectrum Scan

We recommend using this feature only on native SCAN models.

Measuring Absorbance or Transmittance of a sample in the wavelength field, at regular intervals (selectable between 0.1 / 0.2 / 0.5 / 1/2/5 nm)

Multi-wavelengths Fotometry

Measuring Absorbance or Transmittance of a sample at one or more set wavelengths (up to 20).

UV Method for DNA/protein Analysis

There are two preset methods for measuring DNA/protein concentration and evaluation of contamination:

- 1) Method 260/280 nm
- 2) Method 260/230 nm

A method based on the previous ones is available with parameters that can be set by the user.

Energy Scan

Measuring emission energy of the lamps to check the efficiency of the sources.

Wine Chromatic Analysis

Measuring Absorbance at 420/520/620 nm and calculating **Color Intensity** and **Shade** according to OIV-MA-AS2-07B guideline (applicable to red and rosé wines)

2. Installation

The procedure for the correct installation of the Easy UV software is shown below.

2.1. Hardware requirements

- Personal computer with Microsoft Windows® operating system
- Pentium processor or later

- CD-ROM
- No. 2 USB port
- Memory RAM minimum 256MB
- Free space on hard-disk 50MB

2.2. Easy UV software installation

**For computers connected to a company network, access as Administrator may be required.
In this case, contact the network administrator (CED).**

- Put EasyUV disc in the CD-ROM;
- Double Click to open the CD-ROM, and then, double click **Setup.exe** (execute as Administrator) which is under the root directory of CD to start installation (Fig.2.2-1), click **Next**;
- Input user's information (Fig.2.2-2), click **Next**;
- Choose install path (Fig.2.2-3), then click **Next**;
- Be sure that the PC is not connected to the PC (Fig.2.2-4), click **Next**;
- Click **Next** (Fig.2.2-5) to install the driver;
- Accept the License with the chek-box (Fig.2.2-6) and then click **Next**;
- Click **Finish** to complete the installation (Fig.2.2-7).

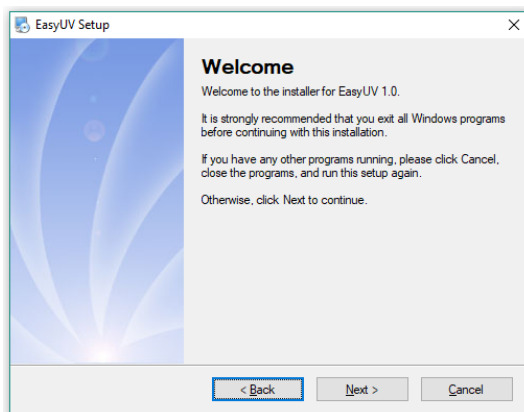
Fig. 2.2-1 Installation box

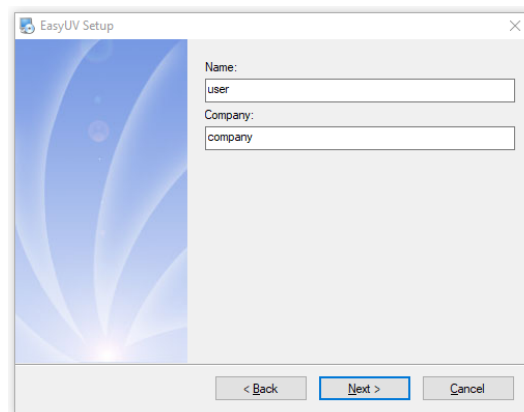
Fig. 2.2-2 User's information

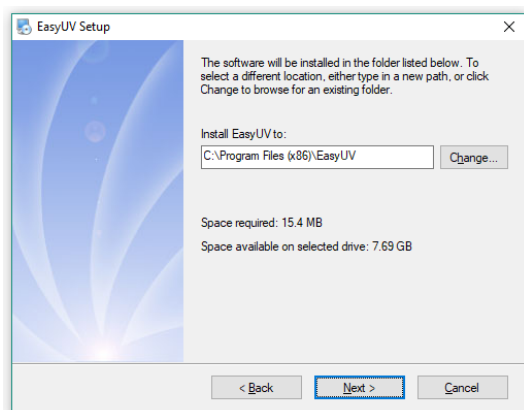
Fig. 2.2-3 Install path

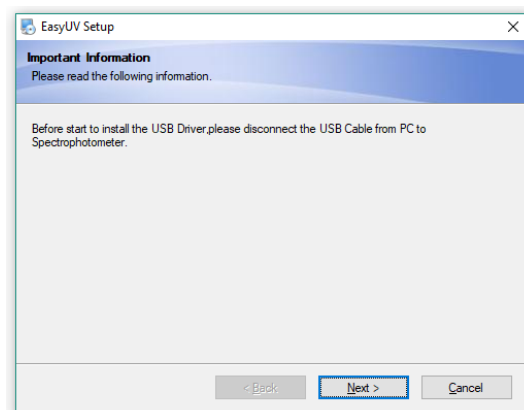
Fig. 2.2-4 Disconnect USB

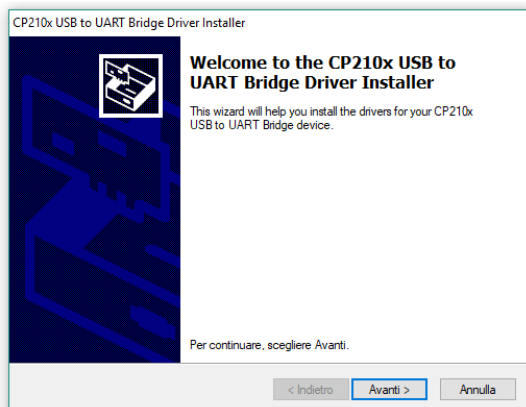
Fig. 2.2-5 Driver installation

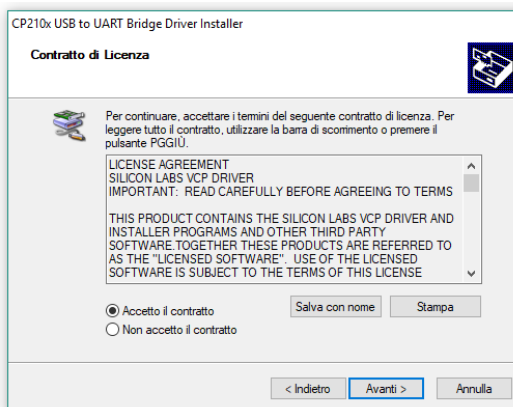
Fig. 2.2-6 License

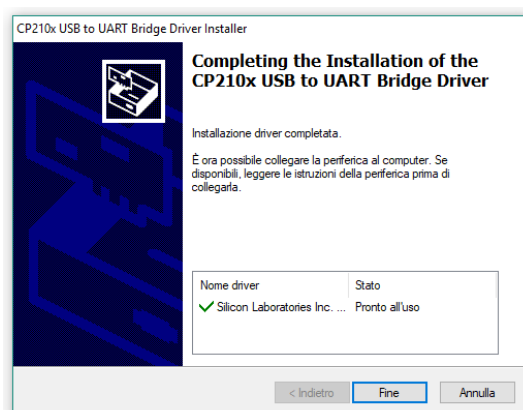
Fig. 2.2-7 Finish

2.3. Spectrophotometer connection

- Connect the spectrophotometer to the PC with the supplied USB cable (Fig.2.3-1) and wait for the automatic configuration of the operating system;
- Click **Finish** (Fig.2.3-2) to exit.

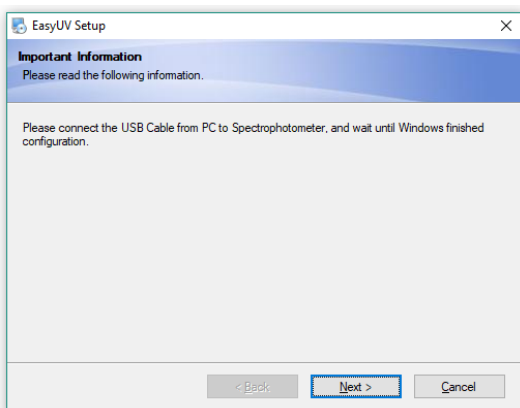
Fig. 2.3-1 Connessione strumento

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2.4. Uninstall

- Menu **Start** → **All file** → **Easy UV** → **Uninstall Easy UV**
- click **Next** to start unistall (Fig. 2.4-1);
- Wait to complete the procedure (Fig.2.4-2), click **Finish**.

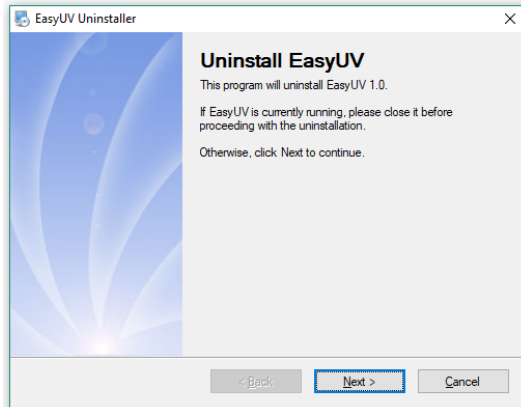
Fig. 2.4-1 Uninstall

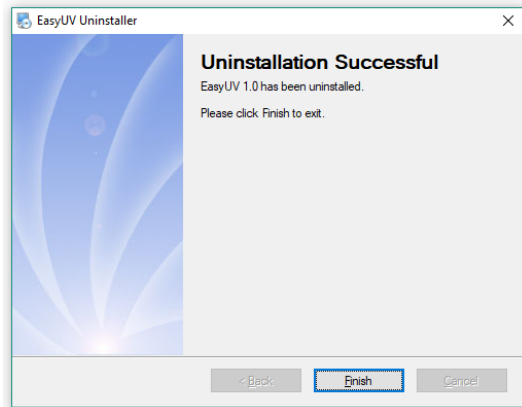
Fig. 2.4-2 Finish

2.5. Check assignment COM port

Be sure the instrument is correctly connected to the PC via the USB cable. Turn on the instrument and wait for the warm-up phase to complete; at the end verify that the display shows the Main Screen HOME (Fig.2.5-1).

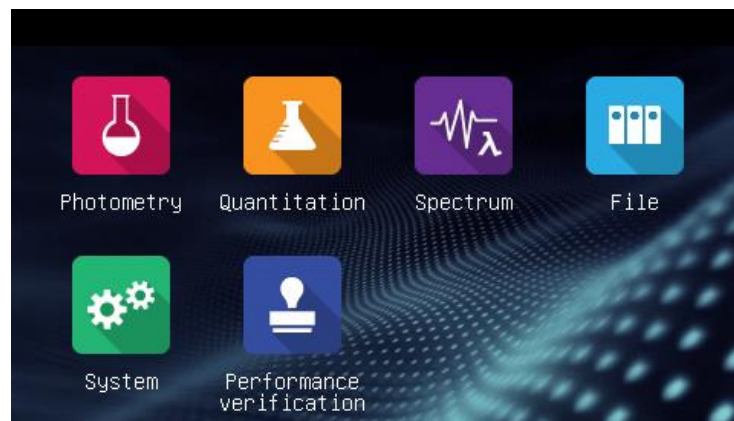
Fig. 2.5-1 Home – Main Screen

From **Control Panel**, go to **Hardware Device**. Open **Ports (COM e LPT)** Fig.2.5-2



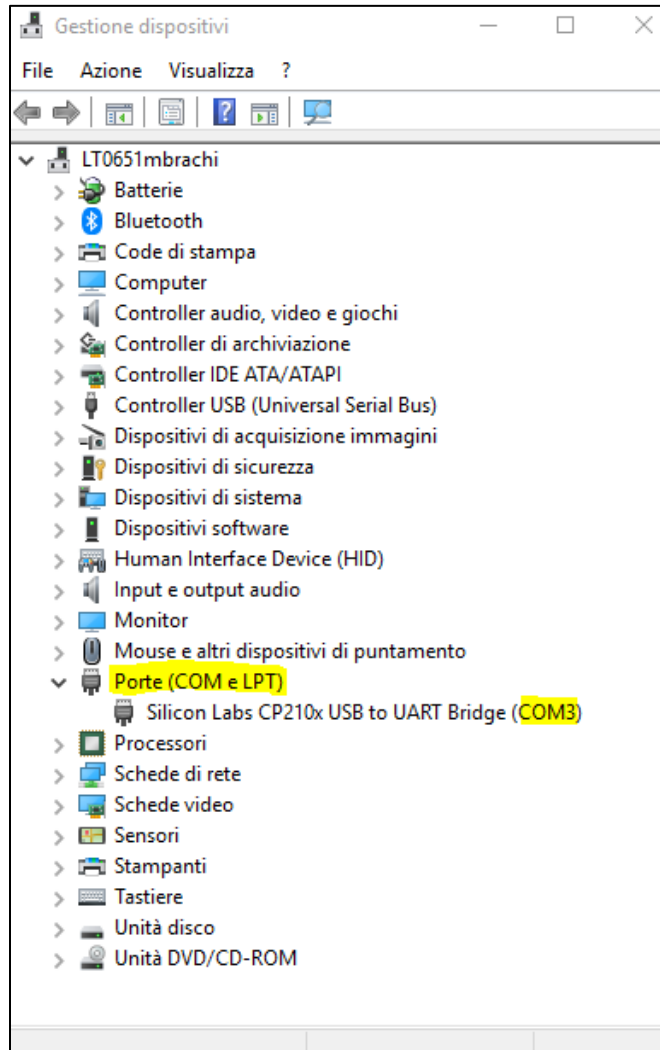
Fig. 2.5-2 Ports COM – Hardware Device Management

Be sure the driver **Silicon Labs CP210x UBS to UART Bridge** is properly installed (no warning signs) and that the assigned **COM Port** is included within the first 8. Otherwise set manually:

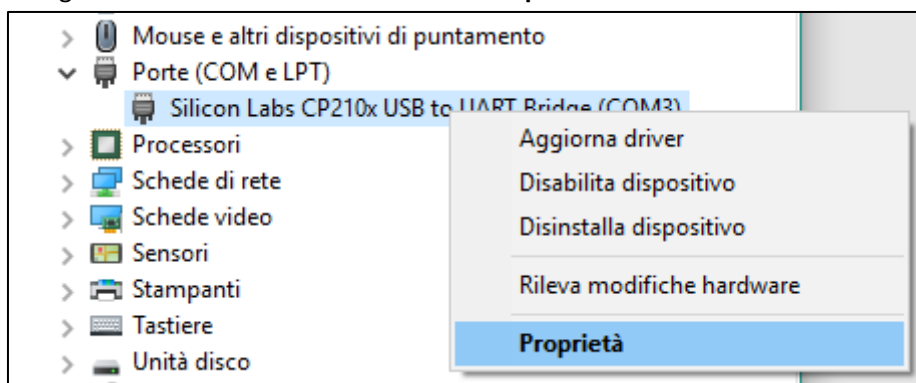
- Right button on **Silicon Labs** and select **Properties**

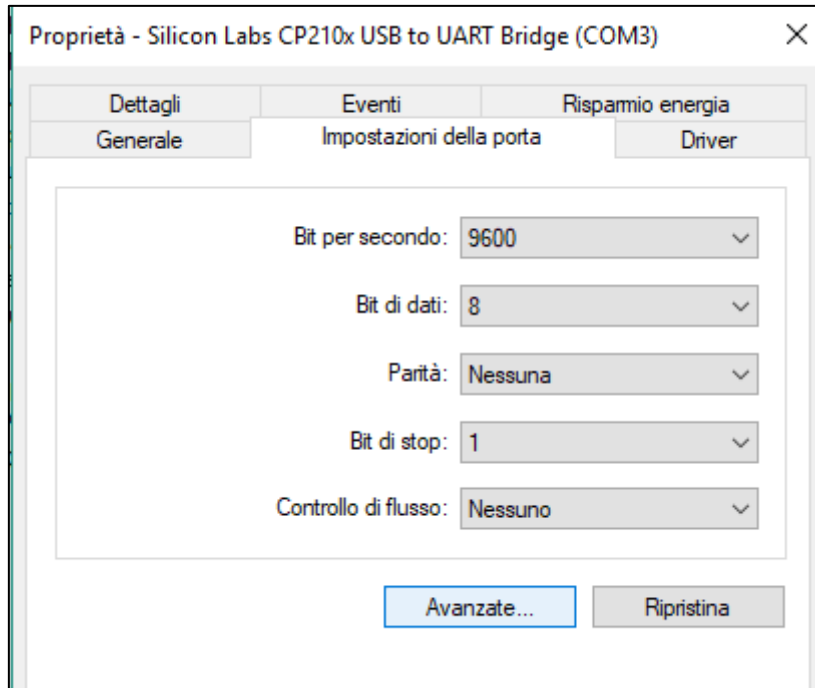- From the **Port Settings** tab click on **Advanced**

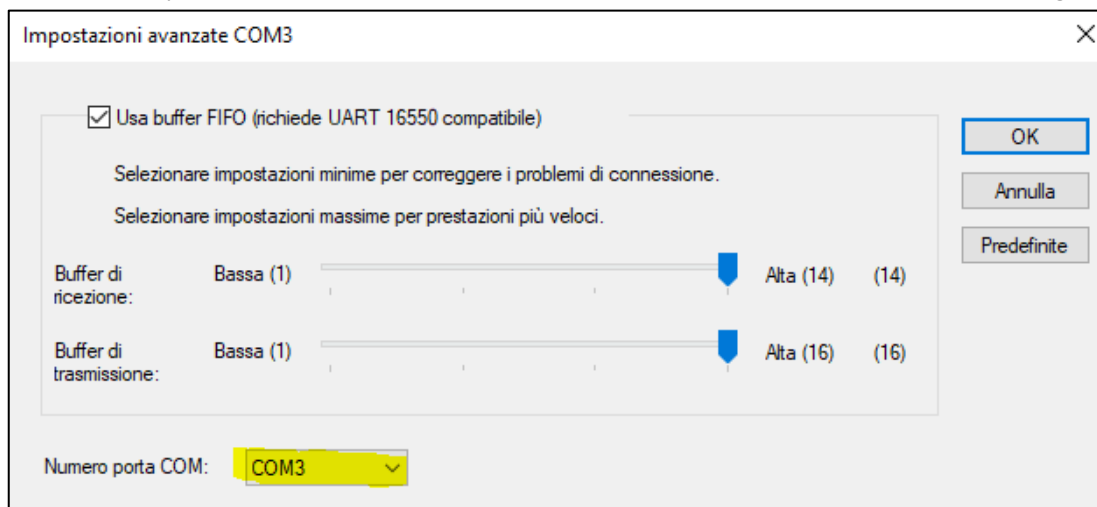- In the **Advanced Settings** tab that opens, in the **COM Port Number** field, select a free COM port between COM1 and COM8. Click **OK** to confirm and exit the Device Manager.

2.6. Starting of Easy UV and first connection

Insert the **UBS Dongle (License for use of the software)** in a USB port on your computer.

Double click on icon  on the desktop to start **Easy UV**.

- menu **View** select **Options** (Fig. 2.6-1);
- **Communication** tab click **Search** (Fig.2.6-2);
- Wait for the automatic search for the tool (Fig.2.6-3)
- At the end of research **Device found** (Fig.2.6-4); click **OK** to confirm.

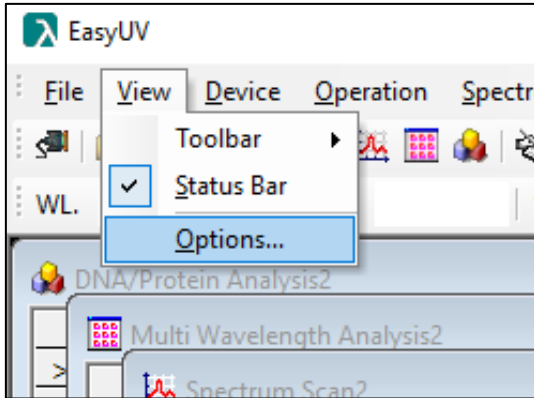
Fig. 2.6-1 Options

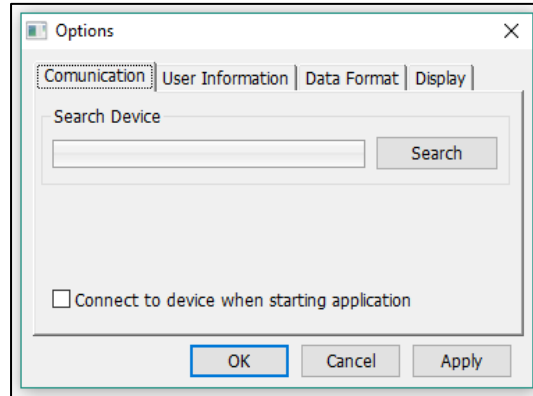
Fig. 2.6-2 Communication tab

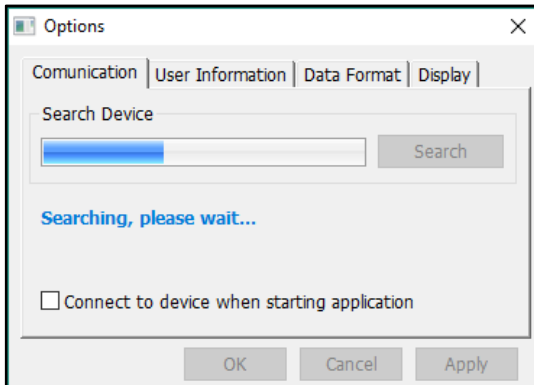
Fig. 2.6-3 Research

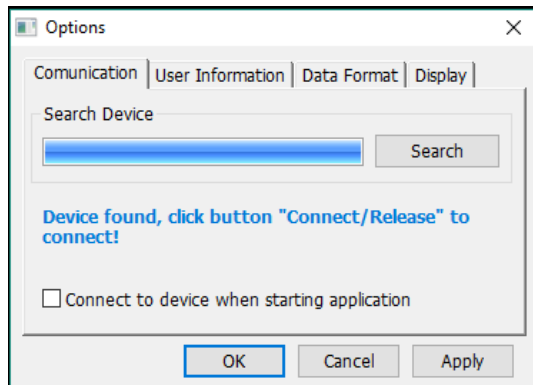
Fig. 2.6-4 Device found

Once the recognition of the instrument has been successfully performed, it is now possible to start the remote control by the program with two modes:

- Menu **Device** → **Connect/Release** (Fig.2.6-5)
- On the Toolbar click on the icon  (Fig.2.6-6)

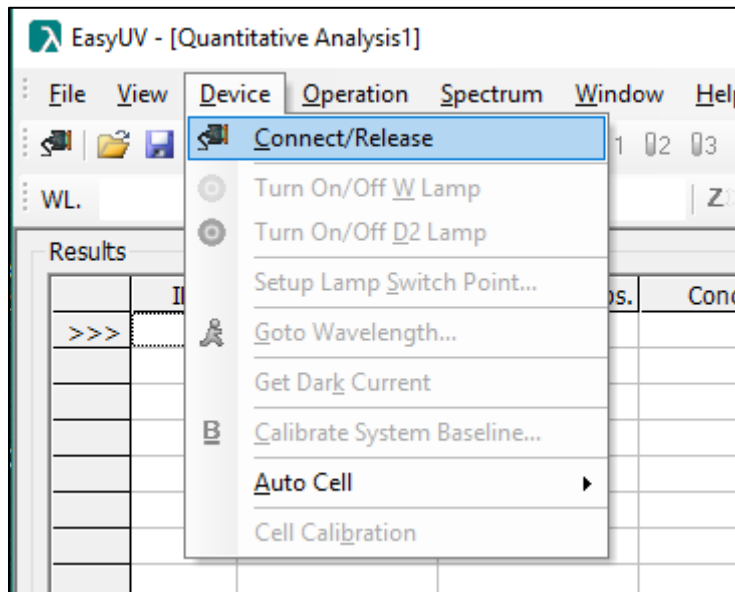
Figure 2.6-5 Connect/Release from Device menu

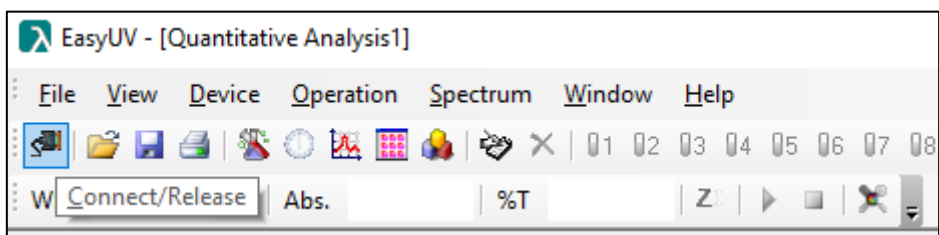
Figure 2.6-6 Toolbar Connect/Release

if the connection has been successful and is operational:

- in the program foot bar the status becomes **Online**
- the function icons are activated in the toolbar
- the spectrophotometer display shows **Online**.

To end the connection click on the icon again .

2.7. User Options Configuration

Menu **View**, select **Options** (Fig.2.6-1). In the box it is possible to configure some settings chosen by the user.

- **User Information tab:** you can select with the check and enter information in the fields that will be added on the printout of the reports (Fig.2.7-1);
- **Data Format tab:** you can choose how the results are displayed (Fig.2.7-2);
- **Display tab:** it is possible to set the software style and the display colors of the screens (Fig.2.7-3);
- Click **OK** to confirm the choices or **Cancel** to esc.

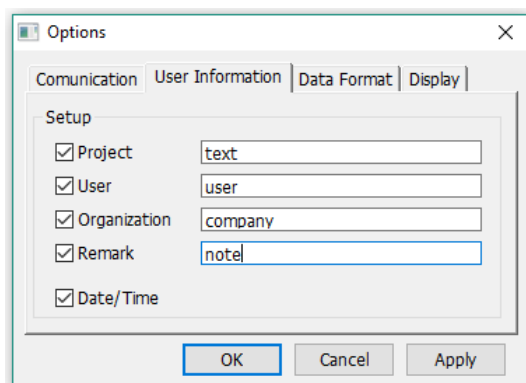
Fig. 2.7-1 User Information

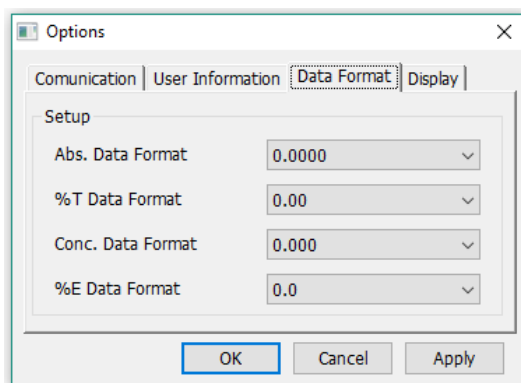
Fig. 2.7-2 Data Format

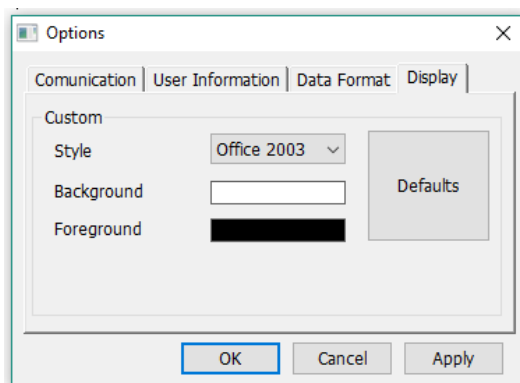
Fig. 2.7-3 Display

3. Introduction

The user interface of the Easy UV software is introduced below.

3.1. Main screen

After starting up Easy UV the main program window is started (Fig.3.1-1), divided into:

- Head bar, which includes the title bar, menu items, and the toolbar;
- Central body (work area), where the windows of the running applications are displayed;
- Foot bar, where the status of the connection is displayed (online / offline), date and time (from the operating system)

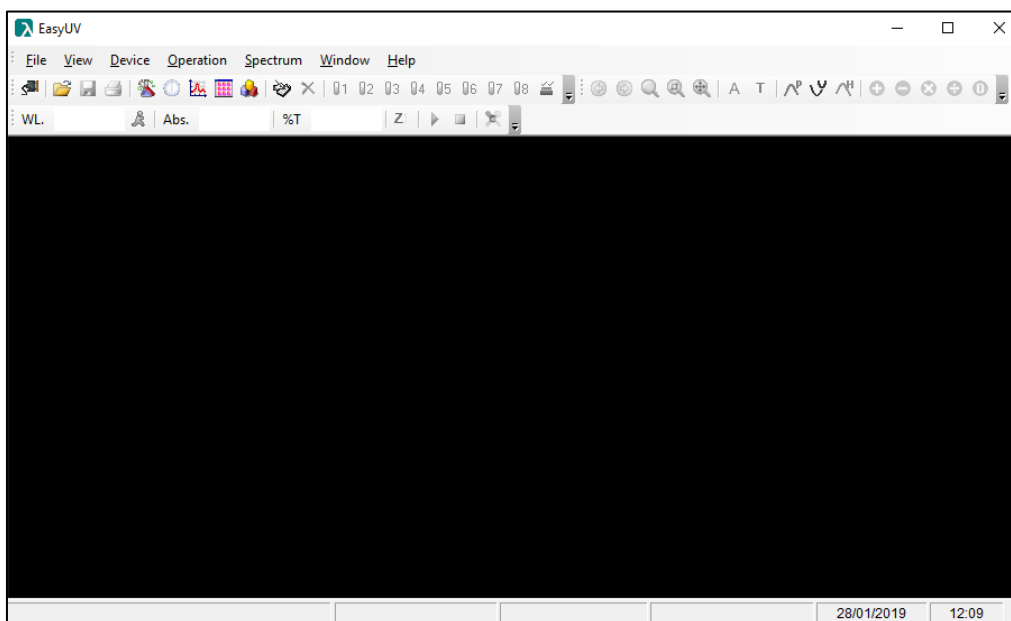
Fig. 3.1-1 Software main screen

3.2. Menu and toolbar

The program offers three modes of use at the discretion of the user:

- Pull-down style Menu (Fig.3.2-1);
- Toolbar (Fig.3.2-2);
- Pop-up menu with right click on the work area (Fig.3.2-3).

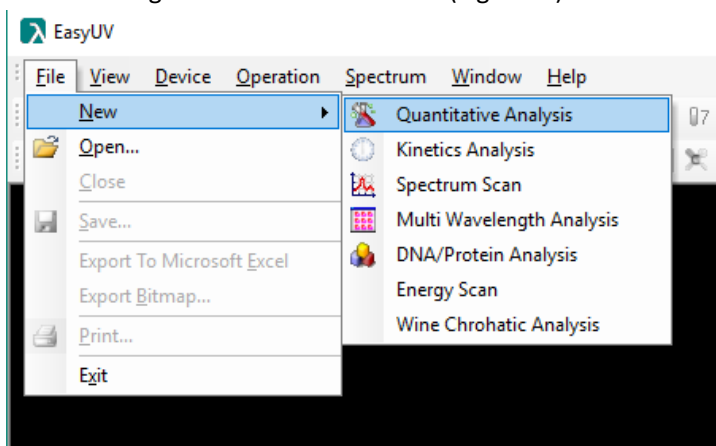
Fig. 3.2-1 Pull-down style menu

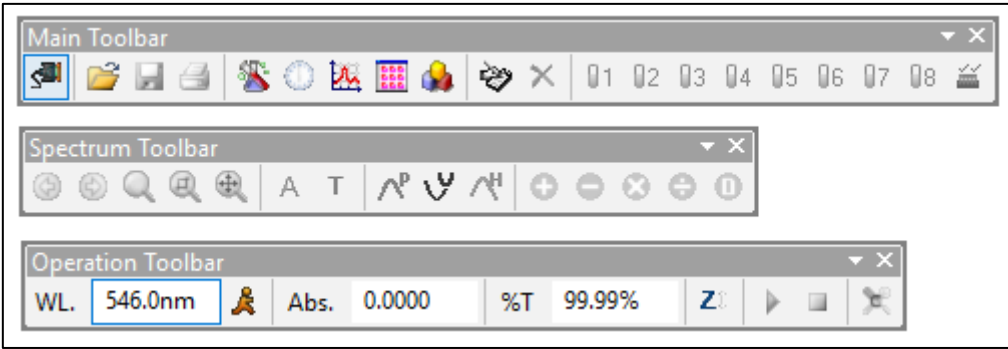
Fig. 3.2-2 Toolbar

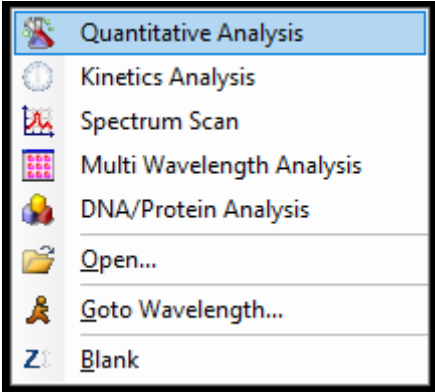
Fig. 3.2-3 Popup menu

3.3. Special functions from menu

EasyUV offers some special features for the use of the user.

3.3.1. Exporting data to Microsoft Excel®

If the Microsoft Excel® software (activated) is installed on your computer, Easy UV offers the user the possibility to export the current window (application) data to an Excel worksheet..

Menu **File** select Export to **Microsoft Excel** (Fig.3.3.1-1).

Wait a few moments and the program creates a spreadsheet in a new Excel file (CartelN) which shows the information related to the test session (*Function and User Information* §2.7) and the measures in the form of a table (Fig.3.3.1-2).

Caution
Depending on the local operating system settings, Microsoft Excel may not automatically recognize the values exported as Number, but consider them Text.
In this case, select the cells containing the values and convert to number format.
You may also need to pay attention to the character used for decimal values (comma "," or dot ".").

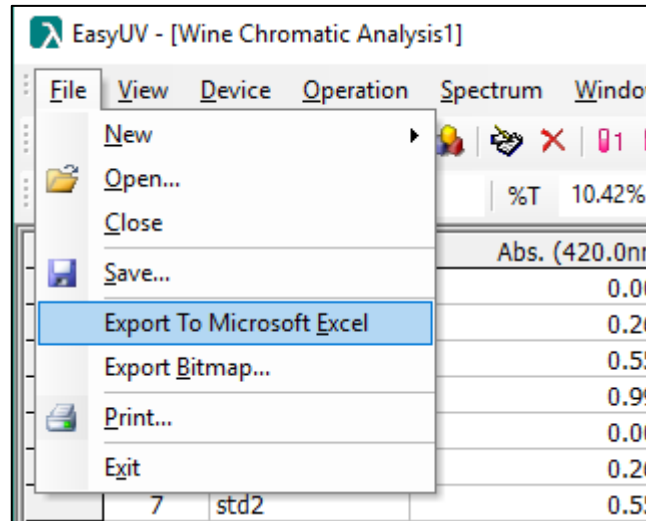
Figura 3.3-1 Export to Excel

 A screenshot of the Microsoft Excel application window titled "Cartel1 - Excel". The spreadsheet contains a report header and a data table. The data table has 7 columns: ID, Sample Name, Abs. (236,0nm), Abs. (257,0nm), Abs. (313,0nm), Abs. (350,0nm), and Abs. (430,0nm). The data rows show three samples, each with an ID and a sample name, and their corresponding absorbance values at the specified wavelengths.

ID	Sample Name	Abs. (236,0nm)	Abs. (257,0nm)	Abs. (313,0nm)	Abs. (350,0nm)	Abs. (430,0nm)
1	Sample1	0.7468	0.8675	0.2917	0.6452	0.9498
2	Sample1	0.7466	0.8677	0.2917	0.6453	0.9495
3	Sample1	0.7468	0.8674	0.2915	0.6448	0.9485

Figura 3.3-2 New cartel exported

3.3.2. Data export in format *.CSV

The program offers the possibility to save data as a data table in format ***.CSV (OpenOffice)**.

Select menu **File → Save**

In the dialog box, select the destination directory, type the file name and select the format

Comma Separated Value file (*.csv) in **Save as**

Click **Save** to confirm (Fig.3.3.2-1).

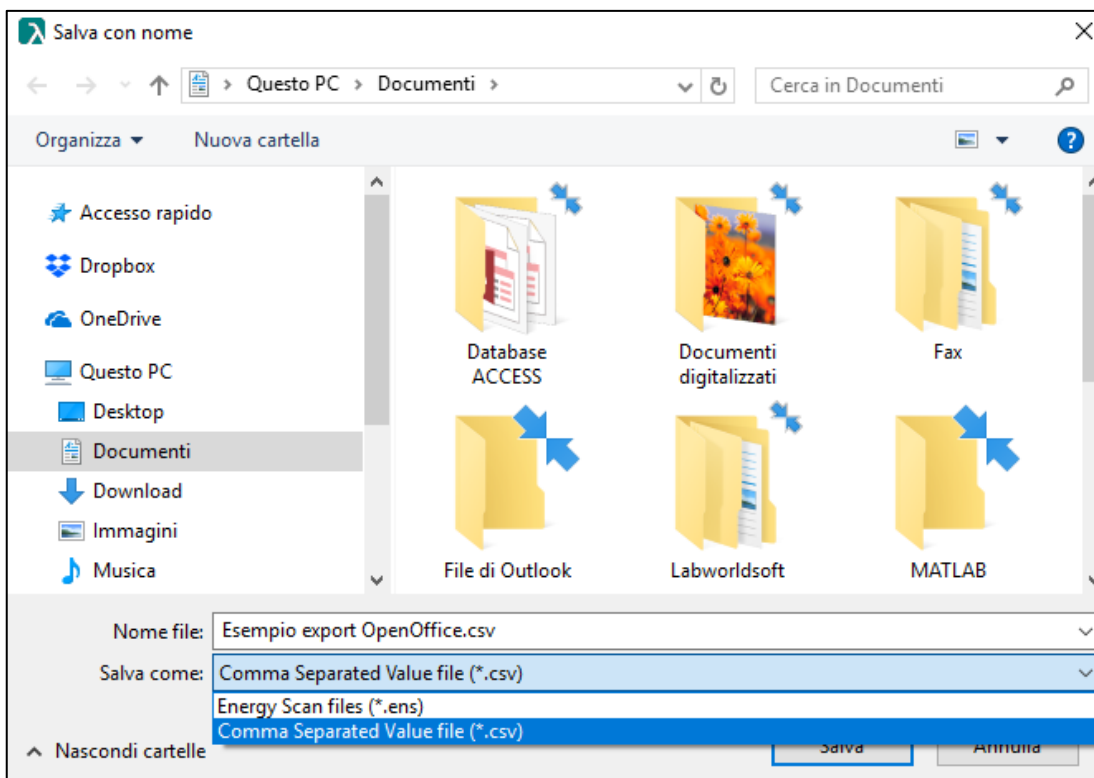
Figura 3.3-3 Saving file *.csv

3.3.3. Menu Device

Menu **Device** offers some setting features of the controlled spectrophotometer. The menu is activated only if the instrument is correctly connected (Fig.3.3.3-1).

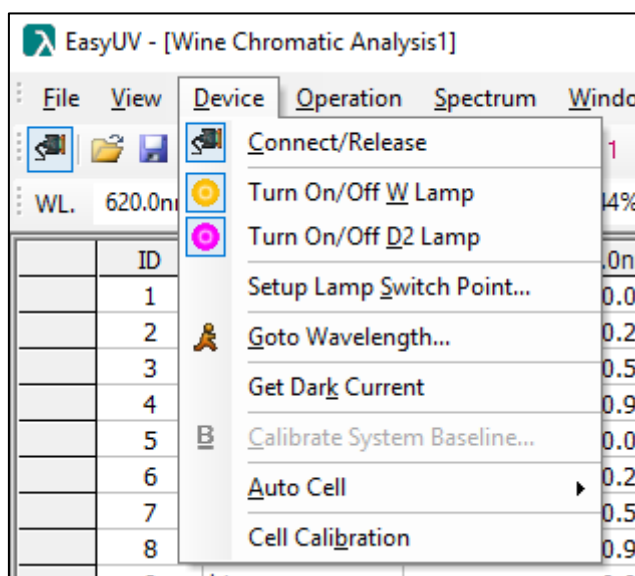
Figura 3.3-4 Menu Device

Connect/Release	Establish connection/disconnection whit the spectrophotometer
Turn On/Off W Lamp	Turn on/turn off the Tungsten lamp (Visible)
Turn On/Off D2 Lamp	Turn on/turn off the Deuterium lamp (UV)
Setup Lamp Switch Point	Setting of the wavelength switch point from Tungsten lamp to Deuterium lamp (Default 340 nm)
Goto Wavelength	Setting of wavelength
Get Dark Current	Perform the measurement of Dark Current

Calibrate System Baseline	Perform the adjustment of System Baseline (only for SCAN models)
Auto Cell	In case of Automatic holder installed on the instrument, Auto Cell manages the positioning of the holder
Cell Calibration	Perform the adjustment of the automatic holder

3.3.4. Functionality with Automatic Holder

In the case of automatic cell holder installed on the spectrophotometer, the software offers the possibility to check the positions of the cell holder.

On the toolbar the menu (Fig.3.3.4-2), the position icons and the setup icon of the cell holder get colored and activated (Fig. 3.3.4-1).

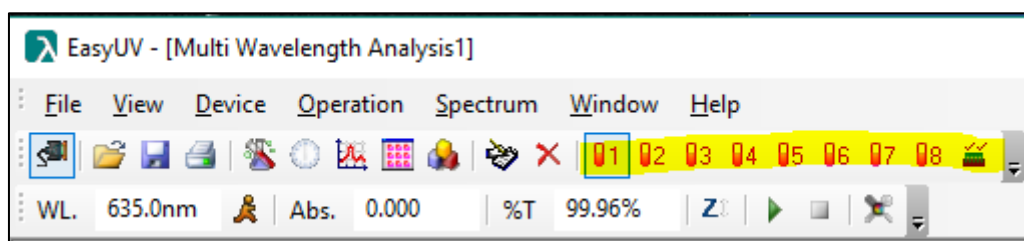
Figura 3.3-5 Automatic holder toolbar icons

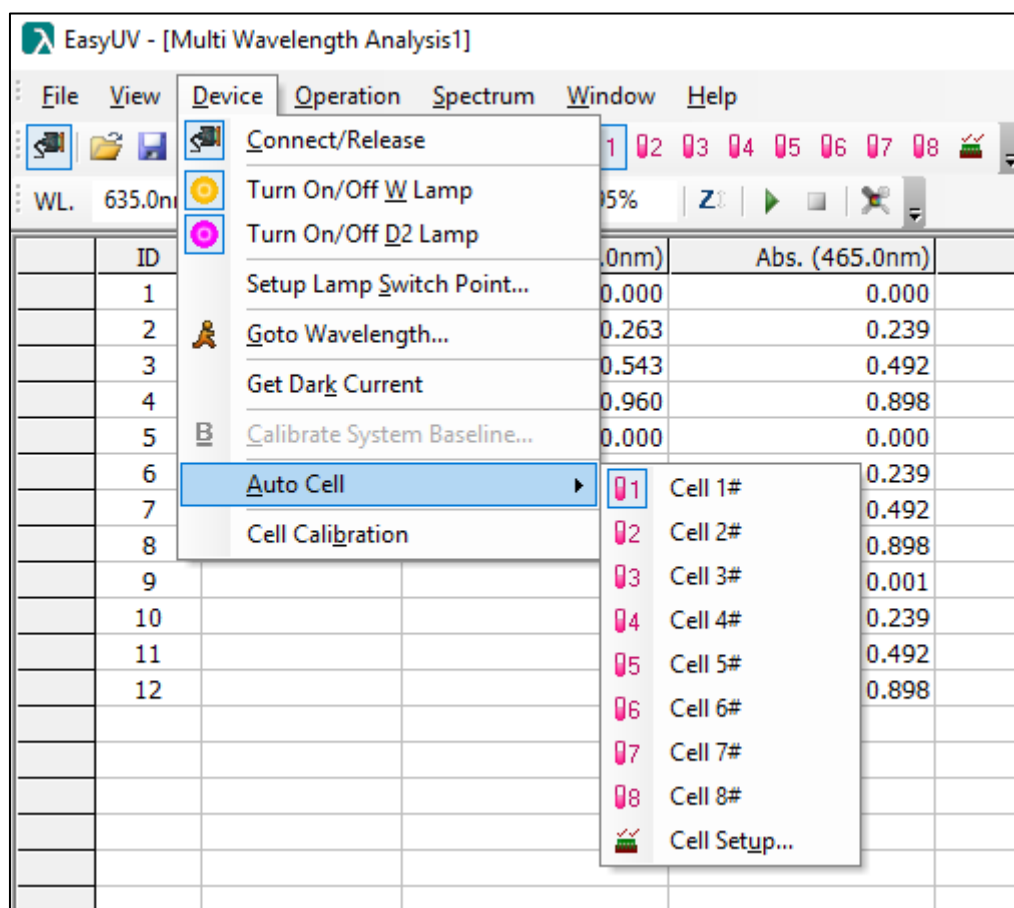
Figura 3.3-6 Automatic holder menu

You can select the positions of the measuring cells by clicking the icon **Cell Setup** (Fig.3.3.4-3); select the cells by ticking.

By checking the **Ref** box, the program takes cell 1 to the measurement position when zero is performed by clicking on the **Zero** icon.

the toolbar or from the pop-up menu.

The worksheet **Quantitative Analysis** consists of two sections (Fig.4.1.-1):

- **Results:** measurement results area
- **Standard Curve:**
 - graphically and analytically displays the calibration curve
Absorbance/Concentration, the calibration points, indicates the correlation coefficient r of the regression;
 - table of the levels used for the creation of the calibration curve, numbered progressively, with indication of the measured photometric value and the concentration level set.

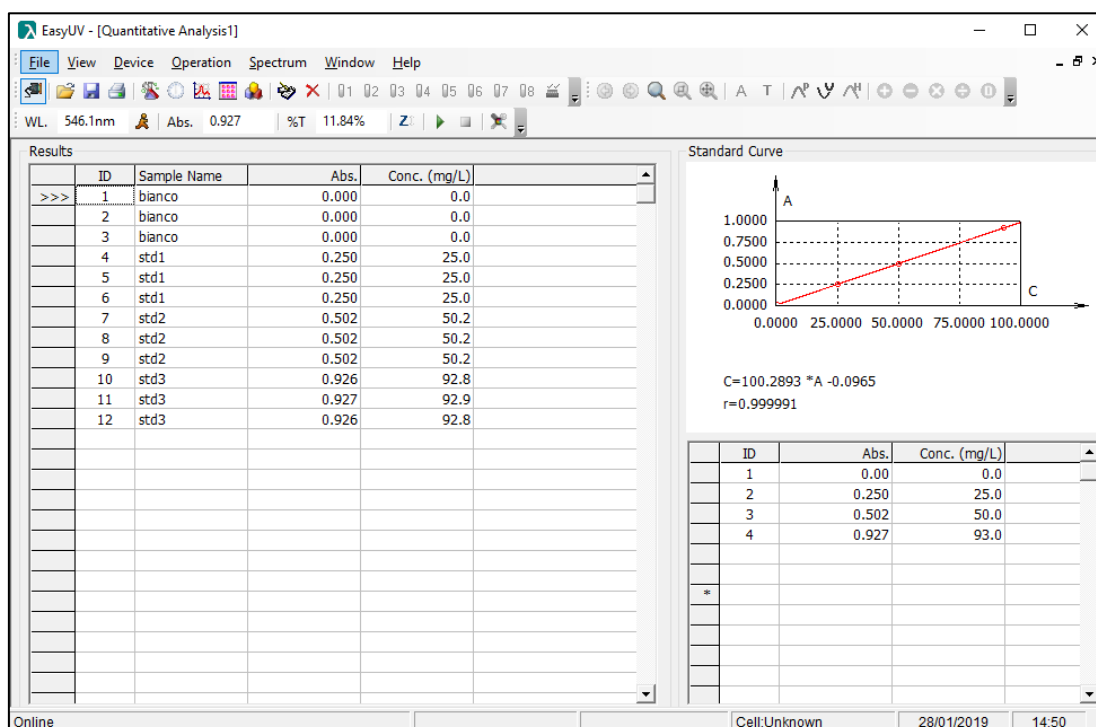
Fig. 4.1-1 Quantitative Analysis

4.1.1. Getting the calibration curve

4.1.1.1. Setting calibration parameters

To access the settings window **Setup** select menu **Operations** → **Setup** or click on the icon .

The Setup window (Fig.4.1.1.1-1) is divided into three sections (identified in columns):

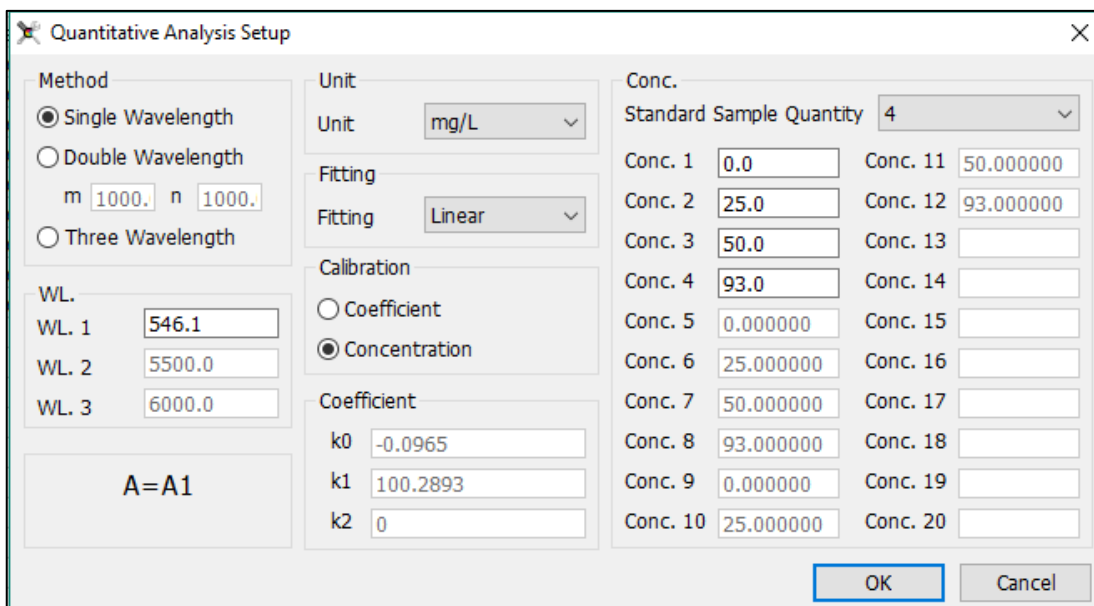
Fig. 4.1-2 Quantitative Analysis Setup

- **Method:** choose the number of desired wavelengths and enter the relative values in the WL fields. The method for calculating the measured photometric value is displayed

Method	Wavelength	Measurement mode photometric value
Single Wavelength	WL.1	$A = A(WL.1)$
Double Wavelength	WL.1 m = coefficiente WL.2 n = coefficiente	$A = m \times A(WL.1) - n \times A(WL.2)$
Three Wavelength	WL.1 WL.2 WL.3	$A = A(WL.1) - A(WL.3) - (WL.1-WL.2) \times (A(WL.2) - A(WL.3))/(WL.2-WL.3)$

- **Unit:** select the unit of measure from those available in the list;
- **Fitting:** select the type of regression of the calibration curve between:
 - **Linear:** linear of the first order
 - **Linear with origin:** linear passing through the origin
 - **Squar:** quadratic of the second order
- **Calibration:** select the method of constructing the calibration curve between:
 - **Coefficient:** allows you to enter known coefficients (factors) without taking standard solution measures; if selected, the editable K fields are activated in which to insert the coefficients in the form $Concentration = f(Absorbance)$:
 - **K0 = intercept (distance from the origin)**
 - **K1 = coefficient of the first degree term (Linear)**
 - **K2 = coefficient of the second degree term (quadratic)**
 - **Concentration:** use of standard solutions with a known concentration to construct the calibration points; if selected, the Conc. column is activated, in which it is possible to choose the number of standard solutions to be used for calibration, from a minimum of 2 to 20 (including any repeated ones). The fields of Conc. In which to insert the relative concentration values with a pre-established order become editable. In this same order the standard measures to build the calibration curve will be carried out.

Click **OK** to confirm the choices made.

4.1.1.2. Calibration with coefficient insertion (Coefficient method)

If the **Calibration: Coefficient** choice was made in the *Setup* and the values of the K coefficients have been entered, the worksheet displays the graph of the calibration curve and the set equation. The standard sector remains empty (Fig.4.1.1.2-1). The Program is ready to perform sample measurements (§ 4.1.2).

The newly created method can be saved to the directory chosen by the user from the menu **File** → **Save**.


Figure 4.1-1 Calibration with Coefficient Method

4.1.1.3. Calibration with measurement of Standard solutions (Concentration method)

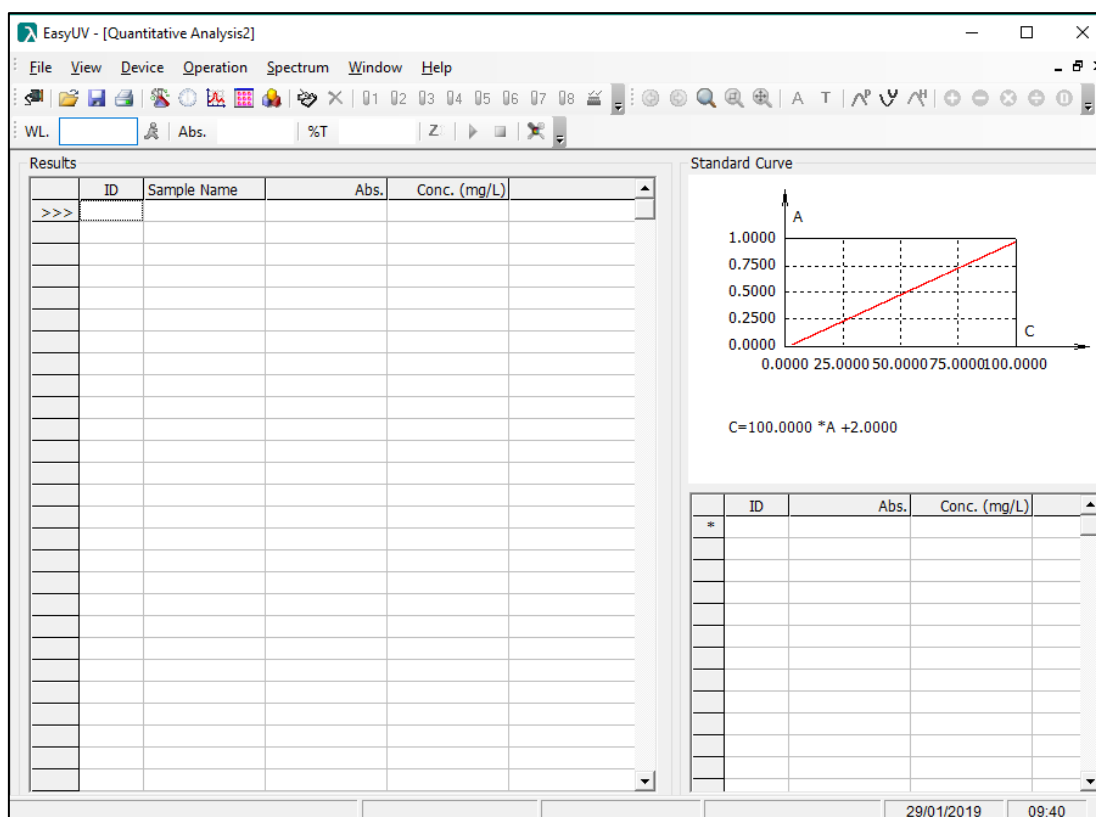If the **Calibration: Concentration** selection was made in the **Setup** and the number of standards has been selected and the concentration values entered, the standard table with the set values is filled in the worksheet (Fig.4.1.1.3-1). To build the calibration curve it is necessary to have the standard solutions of the concentration levels chosen and of the reference (white):

- Place the reference in the cell holder, press the icon **Zero**  ;
- Position the ID1 standard and press the icon **Measure** , the measurement point is displayed on the graph;
- Repeat the previous step for all the set standards.

At the end of the last standard measurement the sheet displays the graph of the calibration curve and the analytical equation of the curve with the correlation coefficient r of the regression (Fig.4.1.1.3-2).

It is recommended to use a statistical approach according to the Eurachem guidelines for the evaluation of the calibration curve and the validation of the measurement methods.

EURACHEM - The Fitness for Purpose of Analytical Methods

possible to measure the samples:

- Prepare the reference in the cell holder and click on the icon **Zero**  ;
- To confirm the correct blanking, click on the icon **Measure**  to measure reference; the result is recorded in the first ID line, the absorbance value of the blank should be as close to 0 as possible. For values greater than 0.01A repeat the Zero and the Measure;
- Prepare the sample in the cell holder and click on the icon **Measure**  ; the data is recorded on the next line.

Repeat the previous point in case of repeated tests and for each sample.

By double-clicking in the **Sample Name** column field, you can type the sample name of the corresponding row (Fig.4.1.2-1).

To save the test session click on the **Save** icon from the toolbar or from the menu **File** → **Save**.

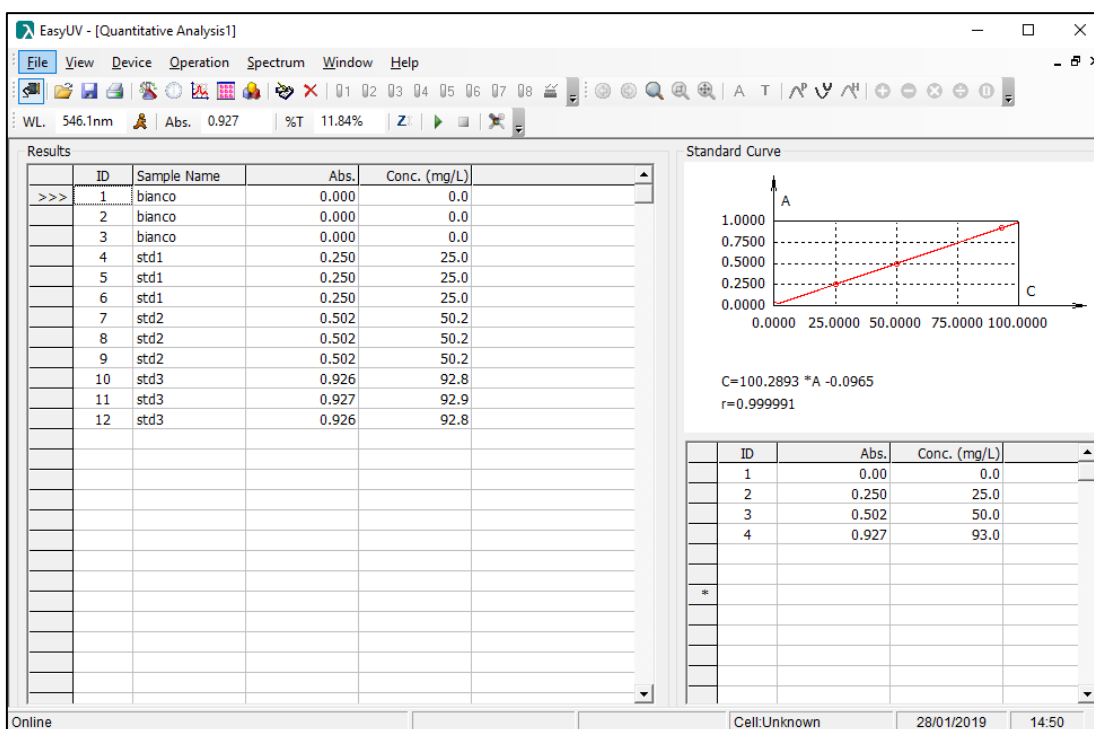
Figure 4.1-4 Quantitative measurement session

4.2. Kinetics Analysis

The Kinetic function measures in photometric mode (Absorbance / Transmittance) a sample over time at regular intervals (selectable between 0.5 and 3600 seconds).

The worksheet displays the Absorbance / time graph and the data table (Fig.4.2-1).

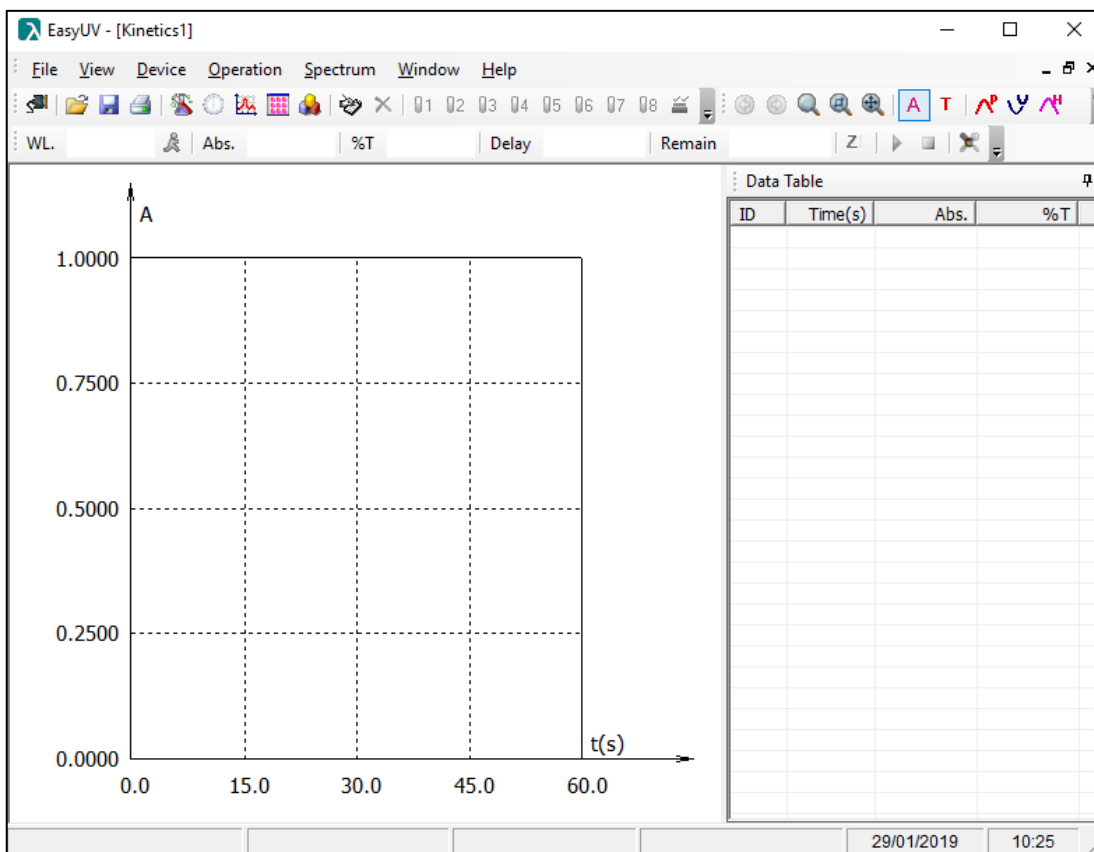
Figura 4.2 Kinetic analysis worksheet

4.2.1. Set measurement parameters

Click on the icon  to access the **Setup** window (see Fig.4.2.1-1). Set the analysis parameters:

- **Photometric Mode:** make the choice to visualize the results between Absorbance (Abs) and Transmittance (% T)
- **Parameter:**

Wavelength (nm)	Enter the measurement wavelength (included in the instrument range)
Delay (s)	Delay expressed in seconds: time between clicking on the icon  and the effective launch of the measures
Total (s)	Time expressed in seconds of the total measurement duration
Interval (s)	Selectable range between 0.5 and 3600 seconds. Time interval in which a measure is recorded

- **Response Mode:** Method of recording the measure *Normal/Slow*
- **Display:** Max = maximum value of the scale of the chart ordinates
Min = minimum value of the scale of the chart ordinates
Check box **Display All Curves:** if selected, in case of several measurements, it shows the curves of the measurements made previously.

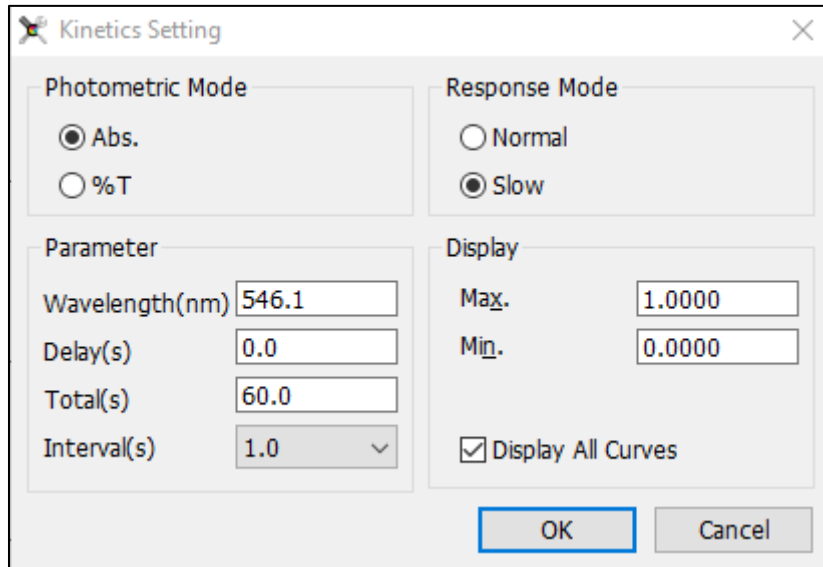
Figure 4.2-1 Kinetic analysis setup

4.2.2. Measure the samples

Once the test parameters have been set up according to the previous paragraph, it is now possible to perform a measurement session.

- Prepare the reference in the cell holder and click on the icon **Zero**  ;
- Prepare the sample in the cell holder and click on the **Measure** icon .

During the measurement, the **Delay** and **Remain** times fields of the toolbar are activated, showing the countdown and the remaining test time respectively (Fig.4.2.2-1).

Figure 4.2-1 Progress Kinetic Analysis

At the end of the measurement, the worksheet shows the graph and the measuring points in the data table (Fig.4.2.2-2)

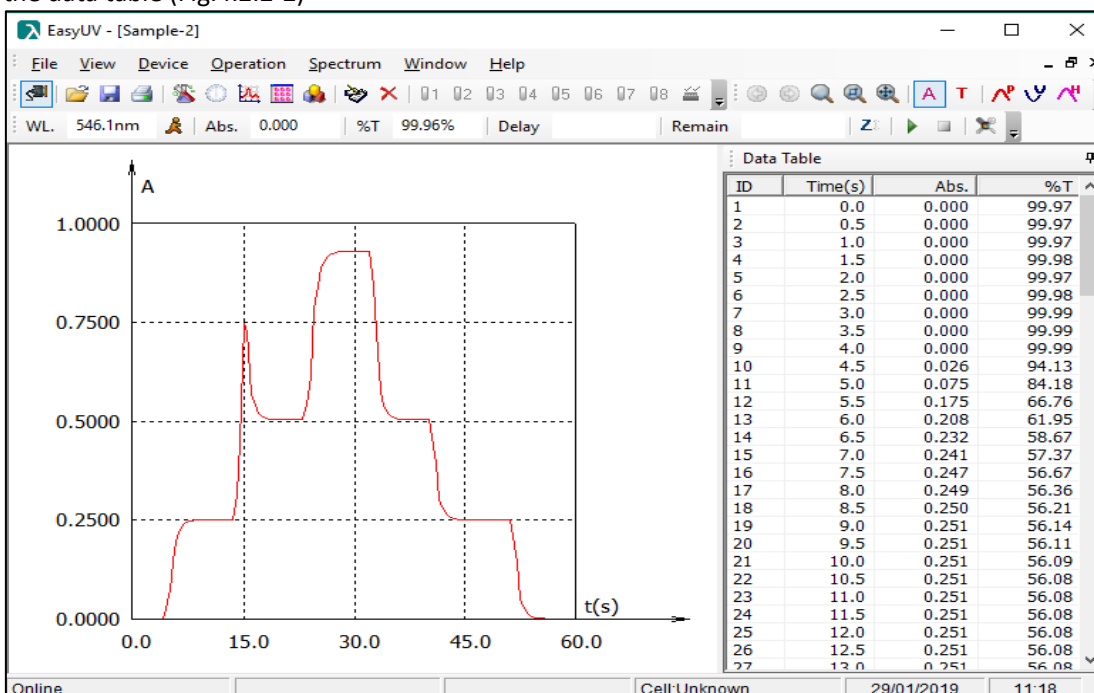
Figure 4.2-2 Kinetic measurement

4.3. Spectrum Scan

We recommend using this feature only on native SCAN models.

The Spectral Scan function measures in photometric mode (Absorbance / Transmittance) a sample in the wavelength field, at regular intervals (selectable between 0.1/0.2/0.5/1/2/5 nm). The worksheet displays the Absorbance / Wavelength graph and the data table (Fig.4.3-1).

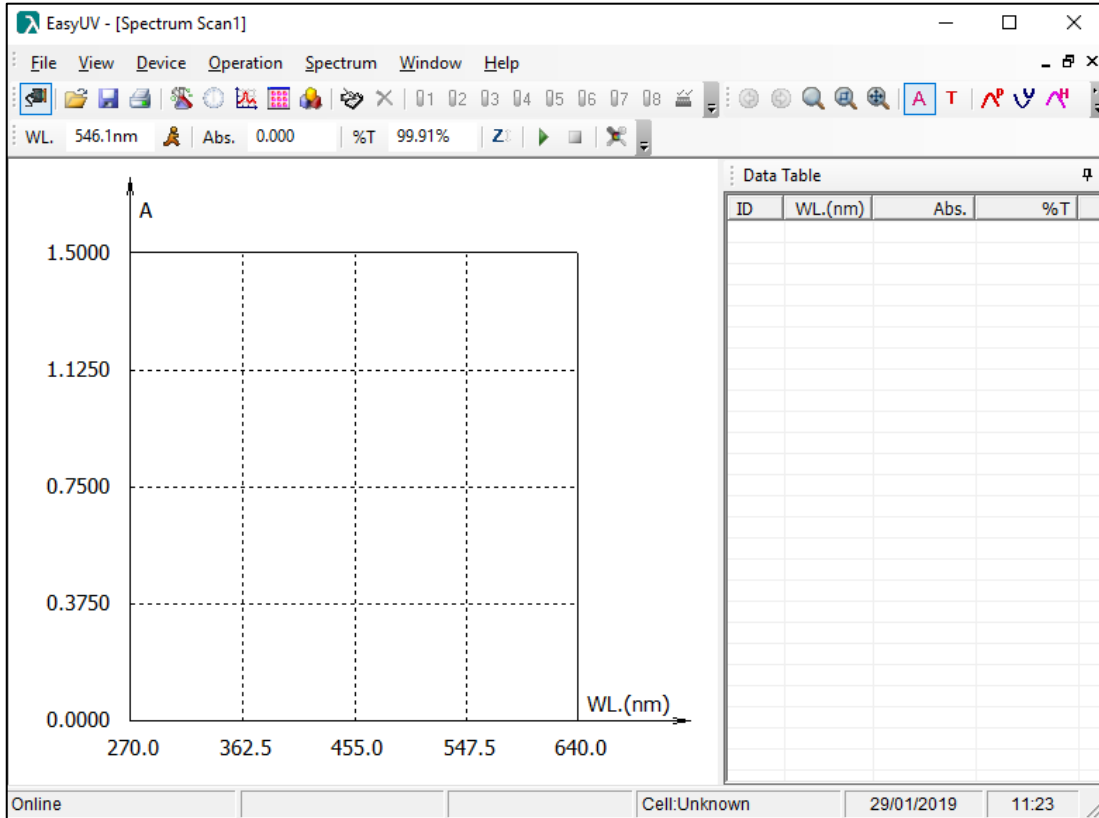
Figura 4.3-1 Spectral Scan Worksheet

4.3.1. Set measurement parameters

Click on the icon  to access the window **Setup** (vedi Fig.4.3.1-1). Set the analysis parameters:

- **Photometric Mode** make the choice to visualize the results between Absorbance (Abs) and Transmittance (% T)
- **Parameter:**

Start (nm)	Enter the initial wavelength (included in the instrument range) higher value
End (nm)	Enter the final wavelength (included in the instrument range) lower value
Interval (nm)	Selectable range 0.1 / 0.2 / 0.5 / 1/2/5 nm. Wavelength range in which a measurement is recorded

- **Scan Mode / Repeat:** Selectable range 1/2/3/5/10 number of automatic repetitions of a measurement
- **Response Mode:** Method of recording the measure *Normal/Slow*
- **Display:** Max = maximum value of the ordinate scale
Min = minimum value of the ordinate scale
Display All Curves check box: if selected, in case of several measurements, shows the curves of the measurements made previously.

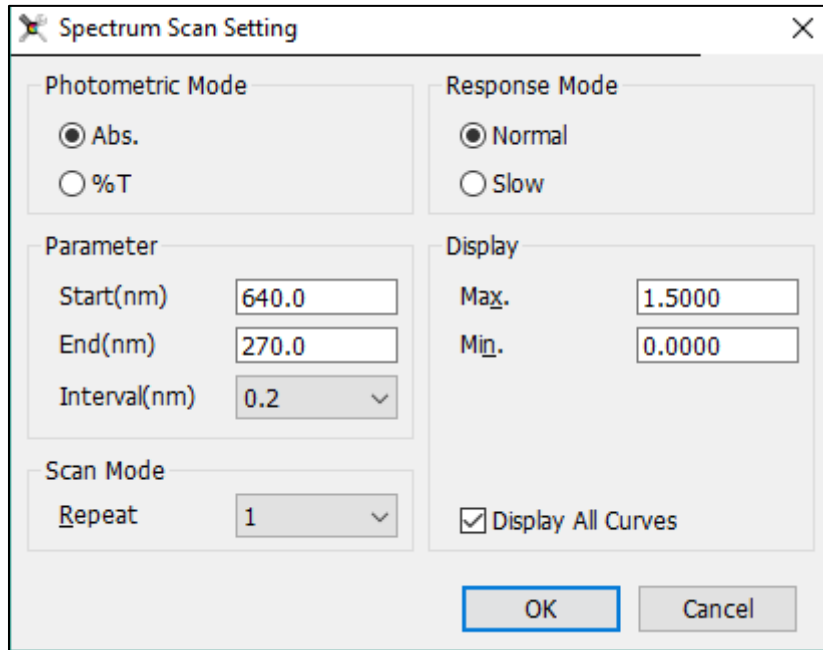
Figura 4.3-1 Spectral Scan Setup

4.3.2. Measure the samples

Once the test parameters have been set up according to the previous paragraph, it is now possible to perform a measurement session.

- Prepare the white in the cell holder and click on the **Zero** icon ; the Baseline window is displayed (Fig.4.3.2-1) with the summary of the measurement parameters, click on Scan to start the construction of the baseline (reference zero) that will be kept in memory for the test session;
- At the end of the Baseline scan, prepare the sample in the cell holder and click on the **Measure** icon .

During the measurement, the **WL** field shows the progress of the wavelengths and the scan curve is displayed in the graph (Fig.4.3.2-2).

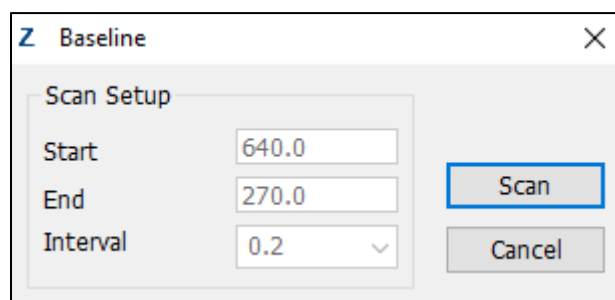
Figura 4.3-2 Confirm Basic Line Parameters (Reference)

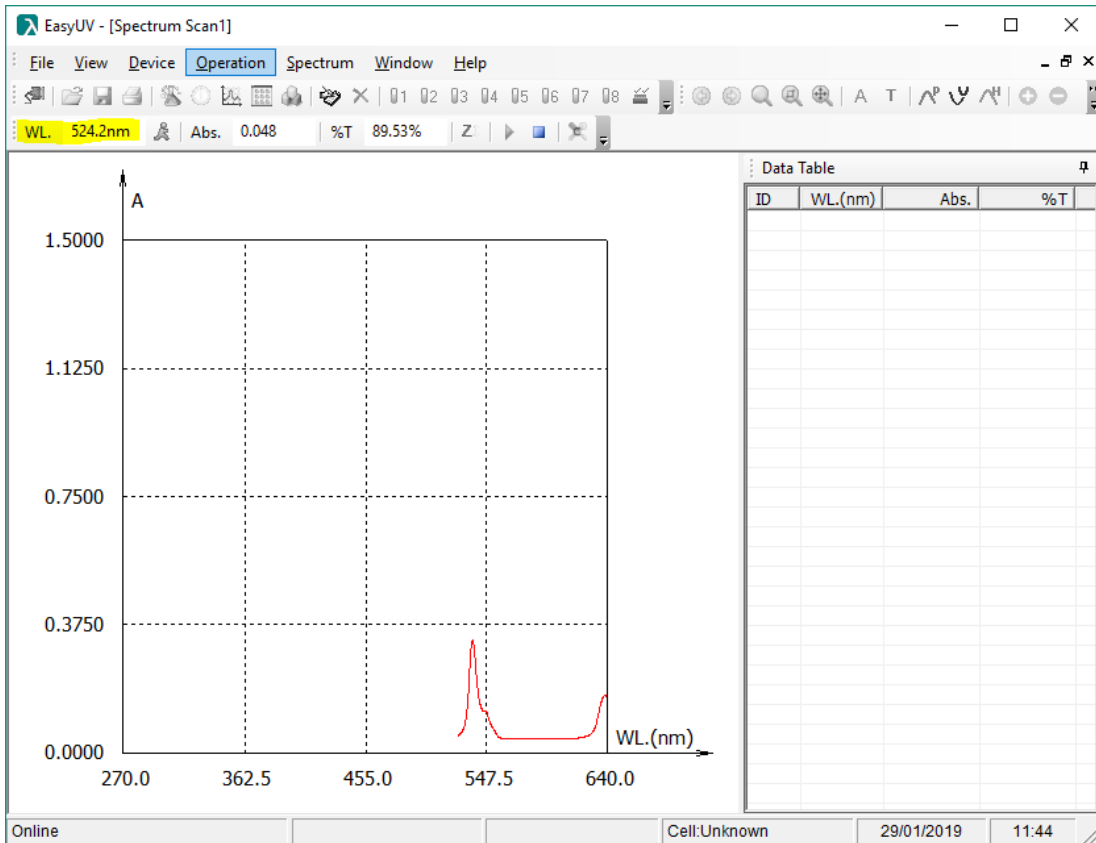
Figura 4.3-3 Spectral Scan Progress

At the end of the measurement, the worksheet shows the complete graph and the measurement points in the data table (Fig.4.3.2-3).

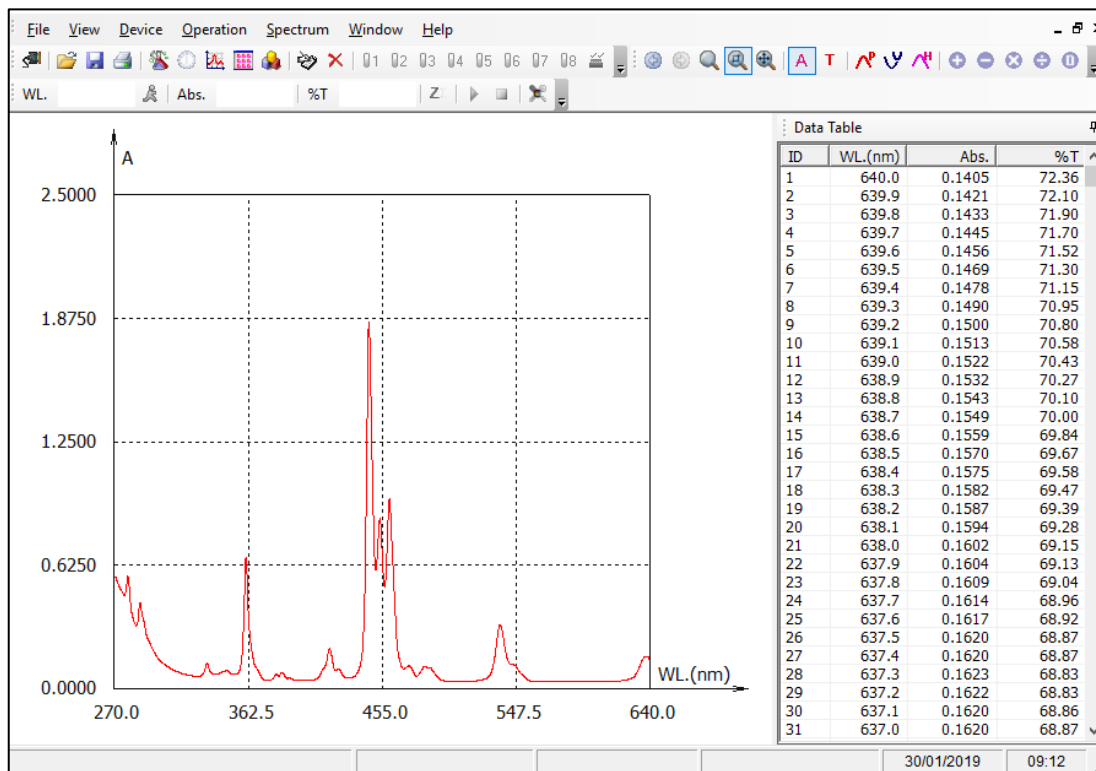
Figura 4.3-4 Spectrum scan

4.3.3. Data processing

At the end of one or more measurements, the program offers some data processing features, accessible from the **Spectrum** menu or from the toolbar (Fig.4.3.3-1).

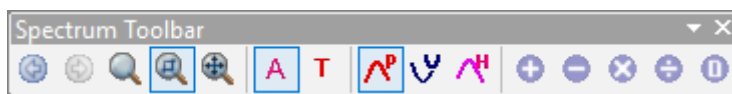
Figura 4.3-5 Spectrum processing tools

Display Mode	Changes the photometric mode display of the Abs /% T data
Current Spectrum	List of Spectrum Scan windows (spectra) loaded
Spectrum Color	List of colors available for displaying the current spectrum
Zoom	set the display of coordinated axes
Peak List	Automatic recognition function of the peaks of the spectrum (Fig. 3.3.3-2). The recognized peaks are numbered progressively and filtered in the data table.
Valley List	Function of automatic recognition of the valleys of the spectrum. The recognized valleys are numbered progressively and filtered in the data table.
Search Condition	Setting the recognition conditions of peaks / valleys. Recommended value between 0.01 and 0.001
Spectrum calculation	Mathematical processing between two spectra (addition, subtraction, multiplication, subtraction) and calculation of the derivative (from first to fourth) of the current spectrum
Smooth	Improve the quality of a spectrum with a lot of noise

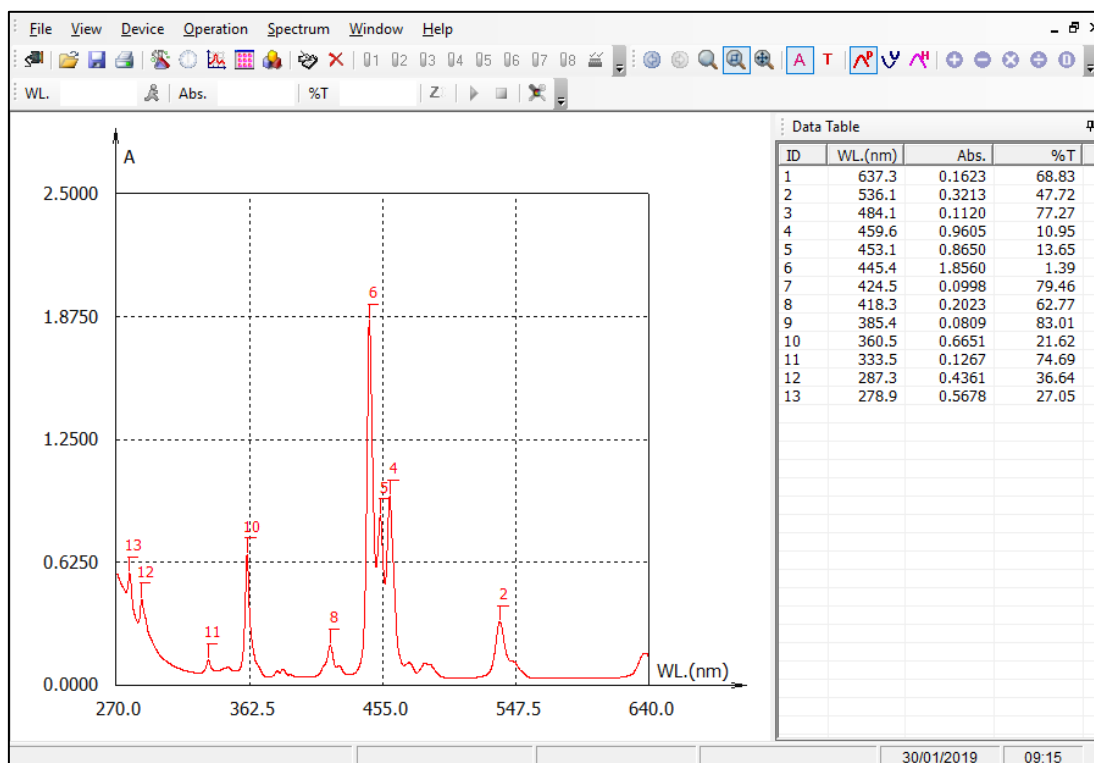
Figura 4.3-6 Display of the peaks

4.4. Multi Wavelength Analysis

The Multi-WL Photometry function measures in photometric mode (Absorbance / Transmittance) a sample at different wavelengths (from one up to 20).

The worksheet displays the measurement results of the samples arranged in a row in a table (Fig.4.4-1).

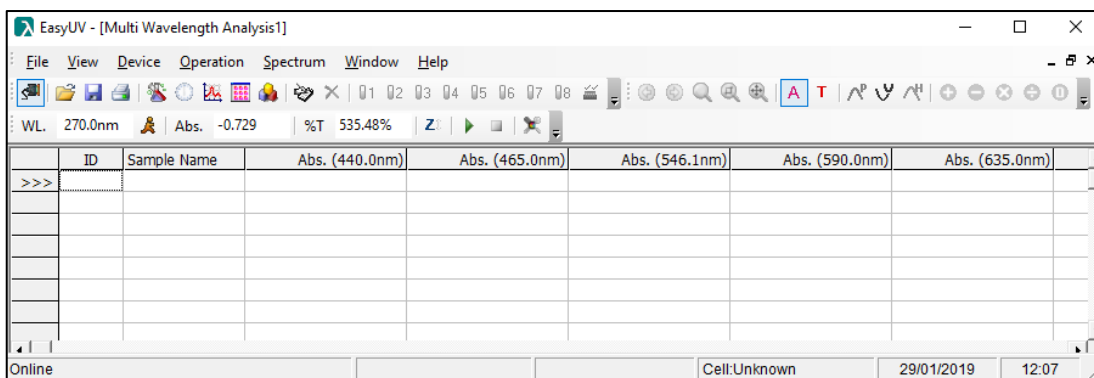
Figura 4.4-1 Multi-WL Photometry Worksheet

4.4.1. Set measurement parameters

Click on the icon  to access the **Setup** window (see Fig.4.4.1-1). Set the analysis parameters:

- **WL No.:** Select the number of wavelengths to measure (from 1 to 20)
- Based on the choice of the number of wavelengths, the relative WL fields are activated in which the measurement wavelengths are to be entered (a maximum of one decimal number is allowed). Press **OK** to confirm.

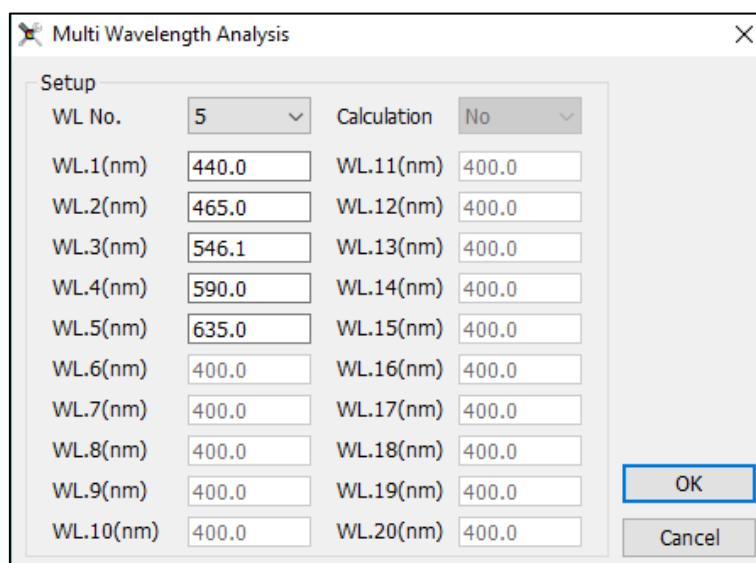
Figura 4.4-1 Multi-WL Photometry Setup

4.4.2. Photometric measurement

At the end of setting the parameters in the previous point, the program is ready to carry out the measurements.

- Prepare the reference in the cell holder and click on the **Zero** icon ;
- To confirm the correct blanking, click on the **Measure** icon  to measure the reference; the result is recorded in the first line, the absorbance value of the blank should be as close to 0 as possible. For values greater than 0.01Abs repeat the Zero and the Measure;
- Prepare the sample in the cell holder and click on the **Measure** icon ; the measured values are recorded by row on the corresponding columns;
- Repeat the previous point in case of repeated tests and for each sample.

By double-clicking in the **Sample Name** column field, you can type the sample name of the

corresponding row.

At the end of the test session the worksheet will be compiled with the tests performed (Fig.4.4.2-1)

To save the test session click on the Save icon from the toolbar or from the menu **File** → **Save**.

ID	Sample Name	Abs. (440.0nm)	Abs. (465.0nm)	Abs. (546.1nm)	Abs. (590.0nm)	Abs. (635.0nm)
1	bianco	0.000	0.000	-0.001	0.002	0.000
2	bianco	-0.001	0.000	0.000	0.001	0.000
3	bianco	-0.001	0.000	0.000	0.001	0.000
4	std1	0.262	0.239	0.250	0.291	0.293
5	std1	0.262	0.239	0.250	0.291	0.293
6	std1	0.261	0.239	0.250	0.291	0.293
7	std2	0.537	0.492	0.502	0.555	0.540
8	std2	0.537	0.492	0.503	0.555	0.540
9	std2	0.537	0.492	0.502	0.555	0.540

Figura 4.4-2 Multi-WL measurement

In case of measures that are considered incorrect, the program offers the user two functions, accessible from the **Operation** menu (Fig.4.4.2-2) or from the toolbar (Fig.4.4.2-3), once the line has been selected of interest:

- **Modify**: repeats the measurement and overwrites the results on the same line;
- **Delete**: Delete the selected row.

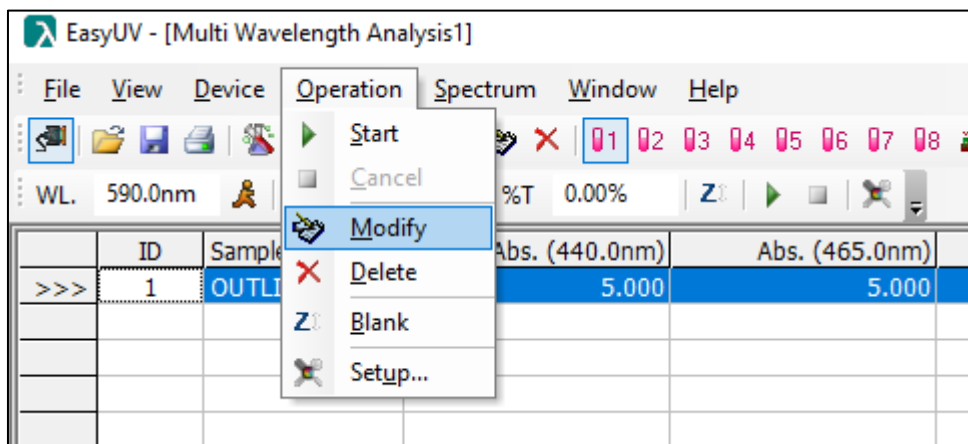
Figura 4.4-3 Edit / delete result

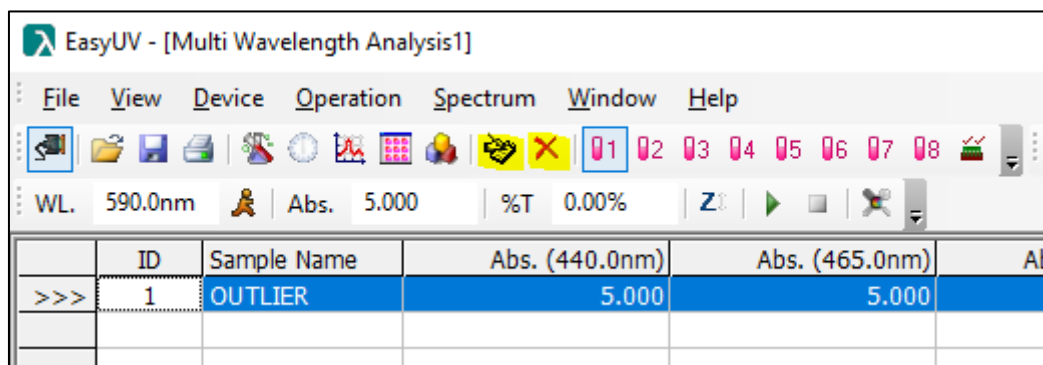
Figura 4.4-4 Edit / delete result icons

4.5. DNA/Protein Analysis – UV Methods

The program offers two preset UV methods for DNA / protein analysis for determining the concentration and / or degree of contamination of the nucleic acid sample from proteins (ratio). The worksheet displays a table whose columns show the absorbance values measured at the set wavelengths, the calculated concentration of nucleic acid (DNA Conc), the concentration of proteins (Protein Conc) and the ratio (ratio) between absorbances measured at the first wavelength and the second (Fig.4.5-1)

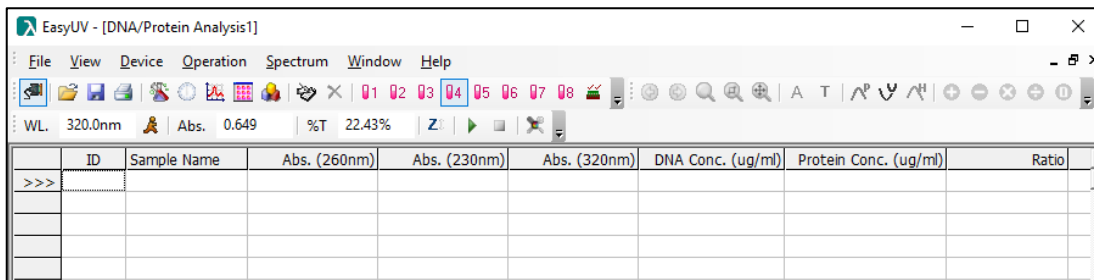
Figura 4.5-1 Worksheet Analysis DNA / Proteins

4.5.1. Set measurement parameters

Click on the icon  to access the **Setup** window (see Fig.4.5.1-1). Set the analysis parameters.

The program calculates the concentrations with the formulas:

- $c_{DNA} = (A_1 - A_{ref}) \times f_1 - (A_2 - A_{ref}) \times f_2$
- $c_{Protein} = (A_2 - A_{ref}) \times f_3 - (A_1 - A_{ref}) \times f_4$
- $Ratio = \frac{(A_1 - A_{ref})}{(A_2 - A_{ref})}$

Method	Wavelengths	Coefficients f _i	Unit of measure
Method 1	WL1 = 260nm WL2 = 280nm Ref = 320 nm (opzionale)	f1= 62.9 f2 = 36.0 f3 = 1552 f4 = 757.3	µg/ml (equivalente a mg/l)
Method 2	WL1 = 260nm WL2 = 230nm Ref = 320 nm (opzionale)	f1= 49.1 f2 = 3.48 f3 = 183 f4 = 75.83	µg/ml (equivalente a mg/l)
User Define	A scelta dell'utente	A scelta dell'utente	A scelta dell'utente

Press **OK** to confirm.

DNA/Protein Analysis Setup

Method

Method 1
 Method 2
 User Define

Wavelength

WL.1(nm) 260.0
 WL.2(nm) 280.0
 Ref.(nm) 320.0

DNA Coefficient

f1 62.900
 f2 36.000

Protein Coefficient

f3 1552.000
 f4 757.300

Conc. Unit

Conc. ug/ml

OK Cancel

Figura 4.5-1 Setup DNA / DNA methods

4.5.2. DNA / protein measurement

At the end of setting the parameters in the previous point, the program is ready to carry out the measurements.

- Prepare the reference in the cell holder and click on the **Zero** icon  ;
- To confirm the correct blanking, click on the **Measure** icon  to measure the reference; the result is recorded in the first line, the absorbance value of the blank should be as close to 0 as possible. For values greater than 0.01Abs repeat the Zero and the Measure;
- Prepare the sample in the cell holder and click on the **Measure** icon  ; the measured values are recorded by row on the corresponding columns, the calculated concentration values are displayed in the Conc fields;

Repeat the previous point in case of repeated tests and for each sample.

By double-clicking in the **Sample Name** column field, you can type the sample name of the corresponding row.

At the end of the test session the worksheet will be compiled with the tests performed (Fig.4.5.2-1)

To save the test session click on the Save icon from the toolbar or from the menu **File** → **Save**.

ID	Sample Name	Abs. (260nm)	Abs. (230nm)	Abs. (320nm)	DNA Conc. (ug/ml)	Protein Conc. (ug/ml)	Ratio
>>> 1	bianco	0.000	0.000	0.000	0.000	0.000	/
2	std1	0.262	0.245	0.102	7.358	14.041	1.119
3	std2	0.861	0.768	0.329	24.593	40.011	1.212
4	std3	1.755	1.535	0.647	51.313	78.518	1.248
5	bianco	0.000	0.000	0.000	0.000	0.000	/
6	std1	0.261	0.245	0.102	7.309	14.117	1.112
7	std2	0.862	0.769	0.329	24.639	40.119	1.211
8	std3	1.758	1.536	0.647	51.456	78.473	1.250
9	bianco	0.000	0.000	0.000	0.000	0.000	/
10	std1	0.262	0.245	0.103	7.313	13.934	1.120
11	std2	0.861	0.769	0.330	24.544	40.087	1.210
12	std3	1.760	1.535	0.647	51.558	78.139	1.253

Online Cell:4 29/01/2019 15:05

Figura 4.5-2 Multi-WL measurement

Wine Chromatic Analysis Setup

WL.

WL.1(nm)

WL.2(nm)

WL.3(nm)

Dilution Factor

Dilution Factor(D.F.)

Color Intensity(I): $I=(A1+A2+A3)*D.F.$

Shade(N): $N=A1/A2$

Figura 4.6-1 Setup Color characteristics of the wine

4.6.2. Measure Color characteristics of the Wine

At the end of setting the parameters in the previous point, the program is ready to carry out the measurements.

- Prepare the reference in the cell holder and click on the **Zero** icon  ;
- To confirm the correct blanking, click on the **Measure** icon  to measure the reference; the result is recorded in the first line, the absorbance value of the blank should be as close to 0 as possible. For values greater than 0.01Abs repeat the Zero and the Measure;
- Prepare the sample in the cell holder and click on the **Measure** icon  ; the measured values are recorded by row on the corresponding columns, the calculated values are displayed in the Color Intensity and Shade fields;

Repeat the previous point in case of repeated tests and for each sample.

By double-clicking in the **Sample Name** column field, you can type the sample name of the corresponding row.

At the end of the test session, the worksheet will be compiled with the tests performed (Fig.4.6.2-1).

To save the test session click on the **Save** icon from the toolbar or from the menu **File** → **Save**.

ID	Sample Name	Abs. (420.0nm)	Abs. (520.0nm)	Abs. (620.0nm)	Color Intensity(I)	Shade(N)
1	bianco	0.000	0.000	0.000	0.0	/
2	std1	0.269	0.263	0.288	0.82	1.022814
3	std2	0.558	0.525	0.539	1.622	1.062857
4	std3	0.990	0.964	0.982	2.936	1.026971
5	bianco	0.000	0.000	0.000	0.0	/
6	std1	0.269	0.263	0.289	0.821	1.022814
7	std2	0.558	0.525	0.539	1.622	1.062857
8	std3	0.990	0.965	0.983	2.938	1.025907
9	bianco	0.000	0.000	0.000	0.0	/
10	std1	0.269	0.263	0.289	0.821	1.022814
11	std2	0.558	0.525	0.539	1.622	1.062857
12	std3	0.989	0.964	0.982	2.935	1.025934

Figura 4.6-2 Measure Color characteristics of the wine

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